

Tec4med

NelumBox  - The worlds most advanced **thermal solution**
for goods requiring temperature control



Operating manual



1	General Information	3
1.1	Identification of manufacturer.....	3
1.2	Scope of delivery.....	4
2	Representation of safety notes	5
3	Notes for the customer	6
3.1	Purpose of the operating manual.....	6
3.2	Liability and warranty coverage.....	6
3.2	Copyright.....	6
4	Basic safety instructions	7
5	Description of the device	8
5.1	Functional description.....	8
5.2	Design and components.....	9
6	Commissioning and operation	10
6.1	Intended use and foreseeable misuse.....	10
6.2	Tips for saving energy.....	11
6.3	General LED / display functions.....	11
6.4	General button functions.....	11
6.5	Connecting / switching on NelumBox®.....	12
6.6	Opening / closing NelumBox®.....	13
	6.6.1 Emergency release.....	14
6.7	Transporting and storing NelumBox®.....	14
6.8	GSM-/GPS-tracking functions.....	15
6.9	Switching temperature profiles.....	16
6.10	Running on power supply.....	18
6.11	Switching batteries.....	19
7	Cleaning and care	21
8	Troubleshooting and disposal	22
9	Technical data	23
10	Declaration of Conformity	24



Thank you for choosing NelumBox®!

In order to ensure a quick commissioning, it is necessary for you to thoroughly read and follow the subsequent notes and descriptions.

1.1 Identification of manufacturer

Name of manufacturer:	Tec4med Lifescience GmbH
Street:	Otto-Hesse-Straße 19 T9
Town:	64293 Darmstadt
Telephone:	+49 [0] 6151 - 360 37 00
E-mail:	info@tec4med.com
Website:	www.tec4med.com



1.2 Scope of delivery



Quantity	Designation
1x	NelumBox®
2x	Li-Ion Battery 
1x	Power Supply 100 - 240 V AC 
1x	RFID card 
1x	Docking-Station 100 - 240 V AC [optional]
1x	Replaceable NIST calibrated Pt-100 temperature sensor
1x	RFID token [optional]

Before putting the unit into operation, check that all parts belonging to the scope of delivery are present.





DANGER!

This symbol signifies **imminent danger** for the **life and health** of persons.



WARNING!

This symbol signifies **potential risks** to **property** and/or the **environment**.



FACTS!

This symbol indicates **important facts** and especially **useful** items of **information**.



Tips for saving energy.



No spray water.



Electronic equipment may not be disposed of with household waste.



Environmentally hazardous.



3.1 Purpose of the operating manual

This operating manual contains important information for safe, correct and economical operation of the device. Observing the instructions will help to avoid dangers, reduce repair and increase the reliability and prolong the service life of the device.

Please keep the manual for future use.

The latest version of this operating manual can be downloaded at any time. Use the following link or QR Code to access the latest version:
www.tec4med.com/nelumboxservice



3.2 Liability and warranty coverage

All information in this manual are provided to the best of our knowledge, on the basis of our experience and findings to date.

The original version of this manual was drawn up in the German language and its technical correctness checked by our company. The translation into the customer's national language/the contractual language has been carried out by a recognised translation agency.

This manual has been drafted with the utmost care. Should you nevertheless establish any gaps and/or errors, please notify us at the abovestated address. Your suggestions for improvements will assist us in designing a user-friendly manual.



3.3 Copyright

All rights are expressly reserved. This manual may not be duplicated or its contents divulged to third parties in any form whatsoever without our written consent.



A basic prerequisite for safe handling and trouble-free operation of this device is acquaintance with the basic safety instructions. It is not permitted to use the device for a purpose other than that stipulated by the manufacturer.

Keep children away from the device at all times!

Never open the housing of the device.

Repair work on the device may only be carried out by Tec4med LifeScience GmbH or by authorised workshops.

Defective parts or components must only be replaced by parts approved of by Tec4med LifeScience GmbH.

Please note the operating and safety instructions of the battery manufacturer RRC.

- If the unit shows visible damage, it must not be put into operation.
- If the connecting cable of this unit is damaged, it must be replaced by the manufacturer or a similarly qualified person to avoid hazards.
- Repairs to this unit may only be carried out by qualified personnel. Improper repairs can pose considerable risks.
- Disconnect the unit from the power supply before each cleaning and care and after each use.
- Medicines may only be stored in original packaging / secondary packaging or suitable containers.
- For infectious material use secondary packaging (approved for infectious or hazardous materials)
- Compare the voltage information on the nameplate with the available power supply.
- Connect the unit only as follows:
with the AC connection cable to the AC network.
- Never pull the plug out of the socket using the connection cable.
- The unit is not suitable for the transport of corrosive or solvent-containing substances.
- The insulation of the unit contains flammable cyclopentane and requires a special disposal process. Dispose the unit properly at the end of its service life.



Danger!



Danger!



Warning!



Danger!



Warning!





5.1 Functional description

The unit is suitable for both mobile and stationary use. It can maintain temperatures below ambient temperature or it can heat goods up. Both uses are viable due to preset temperature ranges.

The unit is a wear-free Peltier cooling [thermoelectric cooling] device with heat dissipation by a fan. The device can be operated at up to 80 % ambient air humidity.





1	upper button	7	cooling/ heating chamber
2	display	8	door
3	cooling grid	9	battery compartment
4	lower button	10	DC connection
5	RFID sensor with green LED light	11	power switch
6	hinged handle		



6 Commissioning and operation

Before commissioning, make sure that the supply cable and the plug are dry.

- If the unit shows visible damage, it must not be put into operation.
- Do not use any electrical equipment inside the unit, unless this electrical equipment is recommended by the manufacturer.
- Do not place the unit near naked flames or other heat sources (heating, strong sunlight, gas oven, heater, etc.).
- Danger of overheating! Always ensure that heat generated during operation can be sufficiently dissipated. Ensure that the unit is at a sufficient distance (at least 10 cm) from walls or objects so that the air can circulate. Do not use the unit in closed systems with a volume of less than 1 m³.
- Make sure that the ventilation openings are not covered.
- Make sure that the loading of the NelumBox[®] does not block the ventilation openings (near the lock & door hinge).
- Do not fill the inner container with liquids or raw ice.
- Never immerse the unit in water.
- Protect the unit and the cables from heat and moisture.
- Avoid throwing or skidding the box or applying excessive external forces.

Operate the device in the following ambient conditions only:

temperature profile:	ambient temperature:
2-8 °C/36-46 °F	min. -10 °C/50 °F; max. 35 °C/95 °F *
15-25 °C/59-77 °F	min. -5 °C/23 °F; max. 45 °C/113 °F *
28 °C/82 °F	min. 10 °C/50 °F; max. 45 °C/113 °F
37 °C/99 °F	min. 17 °C/63 °F; max. 45 °C/113 °F

* Min. 0 °C/32 °F when charging the battery.

6.1 Inteded use and foreseeable misuse

NelumBox[®] is suitable for cooling and heating medical drugs and samples. All materials used are safe for medication. NelumBox[®] may not be checked in as baggage and stored in the cargo hold (Li-Ion batteries). Please check whether the capacity of the unit meets the requirements of the medication you want to keep at the right temperature.





In Australia, Canada, Germany and the United States of America NelumBox® shall not be used for the storage and transport of any type of blood and blood derivatives, organs and tissues.

6.2 Tips for saving energy

A fan inside the unit circulates the air inside the chamber, resulting in better cooling / heating performance.

- Choose a well-ventilated place protected from solar irradiation.
- Do not open the unit more often than necessary.
- Do not leave the unit open longer than necessary.
- Once attached to the power grid, NelumBox® switches to AC-mode. Thus to increase runtime, keep NelumBox® attached to the power grid as long as possible before using the device.



6.3 General LED / display functions

After the selected temperature has been reached, the display turns off and the LED light is illuminated.

The LED light will be on, as long as the display is off and NelumBox® is within the defined temperature range [without error messages].

The display can be activated by pressing one of the buttons. If there is no activity for more than one minute, the display is turned off again (if NelumBox® is still within the defined temperature range).



6.4 General button functions

The buttons can illuminate in the following colors:

- Blue: NelumBox® is cooling, and has not yet reached the target temperature range.
- Red: NelumBox® is heating up, and has not yet reached the target temperature range. / Access via RFID denied. / Battery lower than 5%. / Critical temperature excursion after target temperature was reached.
- Not illuminated: NelumBox® is in the target temperature range.
- Green: Access granted via RFID.



- Yellow: When attempting to open the door via button, but RFID is required. / Battery charge lower than 10%. / Minor temperature excursion after target temperature range was reached. / Door is open for up to 12 seconds.
- White: Door has been open for up to 12 seconds.


6.5 Connecting/ switching on NelumBox®

Make sure that only items or goods that may be stored at the selected temperature range are placed in the device.



Warning!

- Place the unit on a firm base.
- Close the door.
- If the batteries are discharged or the batteries are not used, connect the unit to an AC power socket using the supplied AC power connection cable. Close the door.
- **Note:** The flat part of the plug must point upwards.

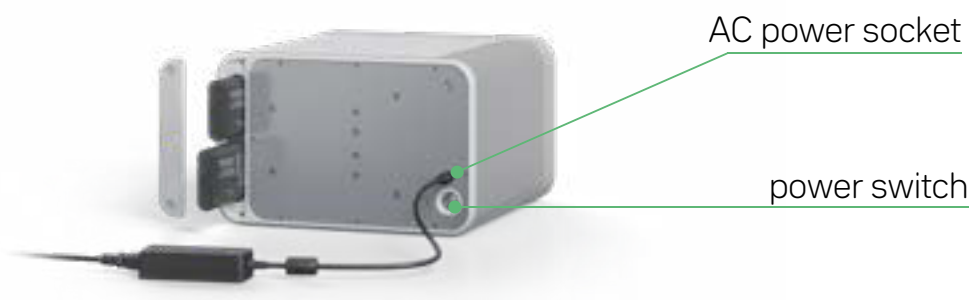

- Press the power switch on the back of NelumBox® to turn on the device. NelumBox® shows the NelumBox® logo during the start-up process.
- The conditioning unit now starts to heat or cool the interior to the previously set temperature range.



After powering up, the battery status and interior temperature of NelumBox® show up on the display. You can also see if NelumBox® was able to successfully connect to the server. Before each transport, make sure that all general conditions for the transport are safely observed!

The batteries can be charged by the Docking-Station. For more information about that device read the RRC2054-2 manual or contact RRC power solutions GmbH.



6.6 Opening / closing NelumBox®



NelumBox® is equipped with an electronic access control using RFID technology [optional]. If the electronic lock is activated, the door can only be unlocked by RFID tokens/cards or via NFC.

NelumBox® can only be opened when switched on [exception: emergency release; see chapter 6.6.1]. To unlock NelumBox®, hold the RFID token/card to the display at the front of the device. Successfully unlocking NelumBox® is confirmed by a green light on the buttons and via the display.

- The lock is unlocked by pulling either the upper or the lower button and by pulling the door. Prevented access is confirmed by the buttons glowing in red and on the display.
- If the electronic lock of NelumBox® is not activated, simply open the device by pulling either the upper or the lower button and by pulling the door.

If unlocking is prevented, do not attempt to open the door by force. This can lead to damage to the device.

- If the unit is opened for more than 12 seconds, a warning signal sounds and the door lights begin to shine in yellow. The signal warns to quickly close NelumBox®'s door to avoid exceeding the set temperature. If the alarm message is ignored, the signal increases in intensity. The display also prompts the user to close the door.
- Close the door by applying pressure to the front of NelumBox®. Press on the area between the display and the hinged handle or on the LED to close the door.



- Do not press on the display and do not push any button while closing the door.
- After successfully closing the door, the audible signal, the lights and the warning on the display disappear.



Warning!



Warning!

6.6.1 Emergency release



In cases of emergency, NelumBox® can also be opened without power supply via cable / battery.

In order to unlock NelumBox® via the emergency release mechanism, push through the sealed opening and trigger the emergency button on the downside of NelumBox® using a suitable device (e.g. a long pen). You can now open the door manually.



6.7 Transporting and storing NelumBox®

- If the unit shows visible damage, it must not be put into operation.
- Danger of overheating! Always ensure that heat generated during operation can be sufficiently dissipated. Ensure that the unit is at a sufficient distance (at least 10 cm) from walls or objects so that the air can circulate. Do not use the unit in closed systems with a volume of less than 1 m³.
- Make sure that the ventilation openings are not covered.
- Protect the unit and the cables from heat and moisture.
- Avoid throwing or skidding the device or applying excessive external forces.
- The permissible maximum overall weight amounts to 15 kg.
- NelumBox® can be put in operation horizontally and vertically.
- To prevent damage make sure the power supply is not plugged in while using NelumBox® vertically.



horizontal operation



vertical operation



- When the GSM and GPS functions are turned off, all other tracking parameters such as humidity, indoor and outdoor temperature, access (time / ID), charge, shock and position are still recorded internally. As soon as NelumBox® reconnects to the Tec4med servers, it automatically sends all retrospectively recorded data to the cloud.

- The same applies if NelumBox® has no connection to the mobile network. In this case NelumBox® shows a missing signal by fading the GPS/GSM symbol.

- As soon as NelumBox® has re-established reception to the mobile network, all retrospectively recorded data is transferred to the cloud.

- When the system is switched on for the first time in a new country, it can take up to 24 hours until the NelumBox® connects to the servers.



6.9 Switching temperature profiles




The control of the unit has a memory function: it remembers the selected temperature range (2-8 °C/36-46 °F; 15-25 °C/59-77 °F; 37 °C/99°F) once the device is shut off.

The previously selected temperature range is set each time the unit is switched on. The selected temperature range is indicated at the top of the display. The basic settings are cold chain „2-8 °C/ 36-46 °F “, ambient „15-25 °C/59-77 °F“ and incubation „37 °C/99 °F“.

Note that it may be intentional if the temperature range cannot be changed. The setting can be changed by authorized personnel or Tec4med.



- You can switch between the different temperature ranges by pressing the upper button for 2 seconds. An acoustic signal and displaying the new temperature range confirms the successfully changed setting. To switch to the next range, it is required to press the upper button for 2 seconds again.

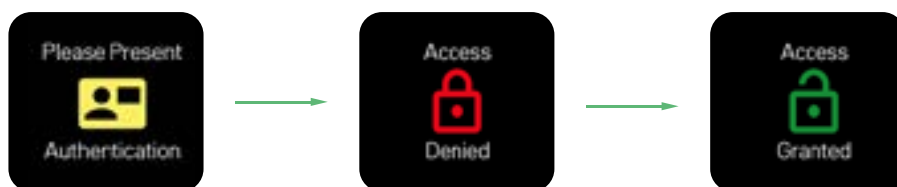

- Access authorization for switching temperature ranges can be granted or denied via the cloud [optional].



- The heating process is indicated by a wave icon.
- The cooling process is indicated by a crystal icon.



- NelumBox® gives a warning (acoustic warning, yellow illuminated door buttons and displaying "Temperature Warning") in case of a minor temperature deviation and an alert (acoustic alert, red illuminated door buttons and displaying "Temperature Alert") in case of a major temperature deviation. Warning and alert can be suppressed by pushing one of the door buttons. Warning and alert will not be given if NelumBox® is approaching the target temperature range for the first time, e.g. after restart or switching the temperature range.



- NelumBox® is operated via the included power supply with the respective country adapter [100 - 240 V AC].



- NelumBox indicates the battery charge level in a battery symbol in the display [white symbol = fully charged battery]. An additional flash symbol is displayed in the battery symbol if AC is connected.
 - If NelumBox® is switched off, it is not possible to charge the batteries. Leave NelumBox® switched on to charge the batteries.
 - The battery status is displayed as a percentage, if NelumBox® is running in battery mode.
-
- An empty battery symbol with a flash symbol is displayed if NelumBox is operated on mains power supply without inserted batteries. Please note that that disconnecting mains power in this state will cause NelumBox® to switch off immediately.
 - For all intermediate states of battery level, the respective battery status is shown on the display. If the charge level is less than 10%, NelumBox® shows a warning in the display, illuminates the buttons in yellow and gives an acoustic warning once per minute. Once the battery charge drops below 5%, NelumBox® shows an alert in the display, illuminates the buttons in red and gives an acoustic alert every 10 seconds. All cooling and heating functions are deactivated, so only the tracking functionalities remain to operate. In this case, please connect NelumBox® to mains power immediately.
-
- The NelumBox®, in the event of a power failure, can guarantee an indoor temperature below 10°C/50 °F for at least 4 hours at an ambient temperature of 25°C/77 °F if two batteries are inserted.



6.11 Switching batteries



To operate NelumBox® properly, the system must either be powered by mains power or by charged batteries. If both batteries are removed at the same time without mains power, NelumBox® immediately switches off completely! Under certain circumstances, this may lead to damage to the loaded goods.

Use NelumBox® only with batteries provided and approved by the manufacturer. The use of other types of batteries can lead to failure or severe damage to the device.



- To change the batteries, open the battery compartment and pull the respective battery out using the tab. [The batteries fit only in one direction.]
- If necessary, turn the lock on the battery compartment lid [e.g. using a coin] to the position where it is possible to remove the lid.



- Batteries can have a different charge level. If two batteries are inserted, the charge level shown on the display is always the average charge level of both batteries.




Warning!



Warning!



- NelumBox® indicates the battery being due for change on the display by indicating either "I" or "II" within the changing icon.
- Please make sure to switch the correct battery.

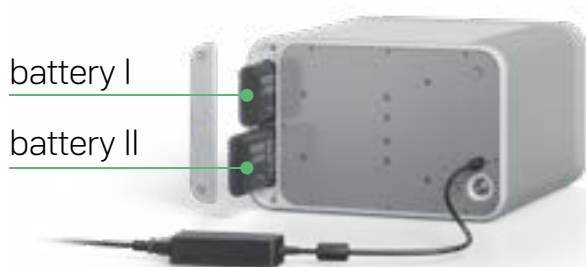

- Do not change the batteries at the same time to maintain the function of the NelumBox®. First replace the first accumulator with a fully charged system before replacing the other accumulator with a fully charged system.
- To avoid failure when replacing the batteries, it is recommended to replace the batteries as soon as they fall below a charge level of 10%.
- Please note, that the batteries must be declared as UN3481 once being packed and delivered by mail. The declaration must state: "Lithium ion batteries in compliance with Section II of PI966".
- Both batteries are fully IATA-compliant and can be carried as hand luggage.





Always disconnect the unit from the power supply or remove the batteries before cleaning and maintenance.

Never clean the unit under running water or in rinse water. Do not use strong cleaning agents or hard objects for cleaning, as these can damage the device.

- Clean the inside and outside of the unit occasionally with a damp cloth.
- Make sure that the ventilation openings of the unit are free of dust and impurities so that the heat generated during operation can be dissipated and the unit is not damaged.



Danger!



Warning!



Electronic equipment may not be disposed of with household waste!

The insulation of the unit contains flammable cyclopentane and requires a special disposal process. Dispose the unit properly at the end of its service life.

The RRC-Li-Ion batteries are part of the GRS-battery system and can be returned at any GRS collecting point globally. Returning the batteries is free of charge! For more information visit: www.grs-batterien.com/

For more information about the web-interface and dashboard visit www.tec4med.com or ask for a manual at help@tec4med.com.

Caution – The battery used in this device may present a risk of fire or chemical burn if mistreated. Do not disassemble, heat above the maximum temperature limit of 45 °C / 113° F, or incinerate. Replace battery with RRC 2054-2 only. Use of other battery may present a risk of fire or explosion.

Dispose of used battery promptly. Keep away from children. Do not disassemble and do not dispose of in fire.

The device and all its accessories must be disposed professionally by the manufacturer. Please also contact the manufacturer in the case of disruptions and emergencies.

Tel.: +49 (0)6151 – 360 37 00

Email: help@tec4med.com

Return to:

TEC4MED LifeScience GmbH, Otto-Hesse-Str. 19, T9,
64293 Darmstadt


If the display shows an error code [E01 - E99] restart the box. If the error code does not disappear or if it returns, contact Tec4med or the source of supply / authorized retailer immediately.



Danger!



Warning!

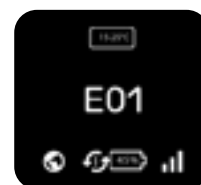



NelumBox



UDI 0 42 6063688 001X XXXXX1

FCC ID:	2AUKO-T4MNLMBXG
Contains FCC ID:	2AJYU-SIM7000G
IC:	25600-T4MNLMBXG
Contains IC:	23761-SIM7000G
HVIN:	T4MNLMBXG
Voltage:	20 VDC 
Power:	100 W
Current:	5 A
Weight:	10,5 kg
W x L x H:	353 x 360 x 225 mm
Net Volume:	4,7 l
IP:	IP20
Ambient temp. charging:	Min. 0°C/32°F; max. 45°C/113°F
Ambient temp. battery mode:	Min. -10°C/14°F; max. 45°C/113°F
Ambient relative humidity:	Max. 80%

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions:
 [1] this device may not cause harmful interference, and [2] this device must accept any interference received,
 including interference that may cause undesired operation.



LABORATORY EQUIPMENT
E505918



Scan code for
service website.



Tec4med LifeScience GmbH
Otto-Hesse-Straße 19
64293 Darmstadt, Germany
+49 6151 360 37 00



NelumBox® is designed and assembled in Germany.



Supplier's Declaration of Conformity

Unique Identifier: NelumBox®

Party issuing Supplier's Declaration of Conformity:

TEC4MED LifeScience GmbH
Otto-Hesse-Str. 19 Block T9
Darmstadt, Hessen 64293 Germany
Telephone: +49[0]6151-360 37 00
Email: info@tec4med.com
Website: www.tec4med.com

Responsible Party U.S. Contact Information / FDA-agent:

Mr. Michael Fritz
Email: fm@bito.com
Ms. Anastasia Wittich
Email: wan@us.bito.com
Company Name: BITO Storage Solutions US, Inc.
Address: 8314 Sherwick Ct/ 20794 Jessup [MD]
Tel.: +1 410 892 6658

Responsible Party Australia Contact Information:

Mr. Bruce Maule
Company Name: Braco Compliance Pty Ltd
Address: Unit 308, 469-481 High St, Northcote, VIC 3070, Australia
ACN/ARBN: 84156023504
Email: admin@bracocompliance.com

Responsible Party Canada Contact Information:

Mr. Ben Hewson
Company Name: Celltech Labs Inc.
Address: 21-364 Lougheed Road, Kelowna, BC V1X 7R8
Tel.: [250] 765-7650 ext. 201
Email: IC-Rep@celltechlabs.com



Innovation, Science and Economic Development Canada ICES-003
Compliance Label: CAN ICES-3 [B]/NMB-3[B]

- This device complies with Industry Canada's license-exempt RSSs. Operation is subject to the following two conditions:
[1] This device may not cause interference; and
[2] This device must accept any interference, including interference that may cause undesired operation of the device.
- Le présent appareil est conforme aux CNR d'Industrie Canada applicables aux appareils radio exempts de licence. L'exploitation est autorisée aux deux conditions suivantes:
1) l'appareil ne doit pas produire de brouillage;
2) l'utilisateur de l'appareil doit accepter tout brouillage radioélectrique subi, même si le brouillage est susceptible d'en compromettre le fonctionnement.
- Antenna Information:
RFID antenna with five windings
Total area: 25x25 mm



NOTE: This equipment has been tested and found to comply with the limits for a Class B digital device, pursuant to part 15 of the FCC Rules. These limits are designed to provide reasonable protection against harmful interference in a residential installation. This equipment generates, uses and can radiate radio frequency energy and, if not installed and used in accordance with the instructions, may cause harmful interference to radio communications. However, there is no guarantee that interference will not occur in a particular installation. If this equipment does cause harmful interference to radio or television reception, which can be determined by turning the equipment off and on, the user is encouraged to try to correct the interference by one or more of the following measures:

- Reorient or relocate the receiving antenna.
- Increase the separation between the equipment and receiver.
- Connect the equipment into an outlet on a circuit different from that to which the receiver is connected.
- Consult the dealer or an experienced radio/TV technician for help.

This device complies with Part 15 of the FCC Rules. Operation is subject to the following two conditions: [1] this device may not cause harmful interference, and [2] this device must accept any interference received, including interference that may cause undesired operation.

The Federal Communications Commission (FCC) warns the users that changes or modifications to the unit not expressly approved by the party responsible for compliance could void the user's authority to operate the equipment.

EU Declaration of Conformity

We,
the company

Tec4med Lifescience GmbH
Otto-Hesse-Straße 19 | T9
64293 Darmstadt
Germany

declare in sole responsibility that the product:

Description: NelumBox®
Type: Gankino
UDI: (01)04260636880018(11)200507(21) (i)

to which this declaration refers, complies with the following directives and standards or other normative documents:

2014/35/EU	Low voltage
2014/30/EU	Electromagnetic compatibility
2014/53/EU	Radio equipment
2011/65/EU	RoHS

EN 61010-1	2010	EN 300 330 V2.1.1	02/2017
EN 61326-1	2013		

DIN 58345	09/2017
-----------	---------

(i): S/N followed by 6 alphanumerical characters uniquely identifying each individual exemplar of the product.

Details of the signatory

Last name: Höler
First name: Nico
Position: Managing Director

Germany
Darmstadt, 14.09.2021
Place and Date


Nico Höler
Signature