

Secure cooling & easy temperature monitoring



Mobile & stationary use




Temperature monitoring



Access control

How to use NelumBox:

1. Press the power switch on the back of NelumBox to turn on the device. The display shows the NelumBox logo during the start-up process. **Cooling starts automatically.**

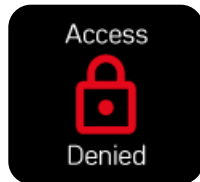

2. NelumBox is equipped with an electronic access control using RFID technology [optional]. The display will show if the access was granted or denied.

3. Open NelumBox by pulling either the upper or the lower button and by pulling the door. Only open NelumBox when access was granted. Do not use any force.



4. Close the door by pressing on the area between the display and the handle or on the door. Do not press on the display. Do not push any button.

Secure cooling & easy temperature monitoring



Mobile & stationary use



Temperature monitoring



Access control

How to use NelumBox:

5. NelumBox constantly cools or heats to maintain the temperature range. The heating process is indicated by the wave icon; cooling by the crystal icon.



6. NelumBox gives a warning (acoustic & yellow display) for a minor temperature deviation and an alert (acoustic & red display) for a major temperature deviation.

7. The battery charge level is indicated in %. An loading flash icon is displayed in the battery symbol if the power supply is connected. **The batteries are now charging.**



power supply

8. The system must be powered by mains power or by at least one charged battery. **Never plug-out both batteries at the same time if no power supply is connected.**

9. NelumBox indicates the battery being due for change on the display by indicating either "I" or "II" within the changing icon. The batteries fit only in one direction.

